Made in the USA

# labSand™

Urine Sample Collection Sand

Noninvasive

100% Natural

Easy to Use

Suitable for most urinalysis tests as a metabolic cage replacement

LabSand is revolutionary biodegradable hydrophobic sand that has a natural non-toxic urine repellent coating. It allows scientists, researchers, and technicians to quickly and easily collect urine samples without resorting to invasive procedures.

#### How to use:

Simply spread the contents of a LabSand bag onto a clean surface and let the animal freely urinate on the sand. The urine will stay on top, allowing you to easily collect the sample with a pipette and vial. Alternatively, you can dip a testing strip directly onto the droplet while it is on the sand. It is that easy!

#### Safety First:

Plastic bags can be dangerous. To avoid suffocation, keep away from babies and children.

**DO NOT EAT LABSAND**

#### EN Urine Sample Collection Sand For Laboratory Use Only

HOW TO USE: Simply spread the contents of a LabSand bag onto a clean surface, and let the animal freely urinate on the sand. The urine will stay on top, allowing you to easily collect the sample with a pipette and vial. Alternatively, one can dip a testing strip directly on to the droplet while it is on the sand. It is that easy!

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## Benefits for Scientists, Researchers, & Technicians:

Easy collection and time saving

100% natural and non-toxic

Noninvasive

Recommended by top pharmaceutical companies

Obtain a clean, uncontaminated sample

Maintains high standards of animal welfare

Biodegradable & easy disposal

Scan to learn more about Lab Sand. Visit [www.labsand.com](http://www.labsand.com) today to learn more about this unique product.

Manufactured by: Coastline Global, Inc. 827 Lincoln Ave, Suite B11,  
West Chester, PA 19380 Tel: 888-999-3050 | [info@coastlineglobal.com](mailto:info@coastlineglobal.com)

Made in the USA